QUICK GUIDE

A/Software installation

Use the "setup" application on the IviumSoft installation CD and follow instructions. IviumSoft will be installed on your C- or D-drive (default) in the "IviumStat" directory. This directory will include example data files, drivers, Labview example and all manuals. Also a shortcut to




2) Pre-Install Drivers:

Navigate to the newly created directory: ..\IviumStat\iviumdrivers and run "Ivium_driverinstaller.exe" to install 2 Ivium instrument drivers. When prompted click "next"; then at the Windows Security pop-up click "Install this driver software anyway"; upon completion click "Finish".



B/Connecting your instrument

- All Ivium potentiostats have a unique serial number that is connected with the device driver. The first time the potentiostat is connected to a specific computer, the driver will be loaded. Normally this will go automatically and take a few moments. If prompted (Windows XP), choose "automatic" installation and to "continue anyway". After the driver has been loaded, the IviumSoft can be started and the device serial number is visible in the readout window at the top left of the user interface.
- The pocketSTAT is USB powered: connect the delivered USB cable between pocketSTAT and PC/Laptop and the instrument is ready to operate.

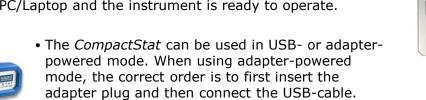

• For all other Ivium-potentiostats, first insert the power cable, then connect the USB-cable. Next, switch the instrument on.



Options - Setting the device and data handling options, incl. FRA options

- HTML help file: content sensitive and topic search

About - Shows IviumSoft version and corresponding .dll [no#]

- Data and method file management.

- Device maintenance and special operations





Close

Scan Chan Olay

Hide all

Show all

☐ Hide previous

All extradata

✓ scan 1:[cycle 1]

Delete

Delete all


Data appear

no labels

with the Ivium designed Electrochemical.


E-Mail: info@ivium.r

- 3. Advanced parameters

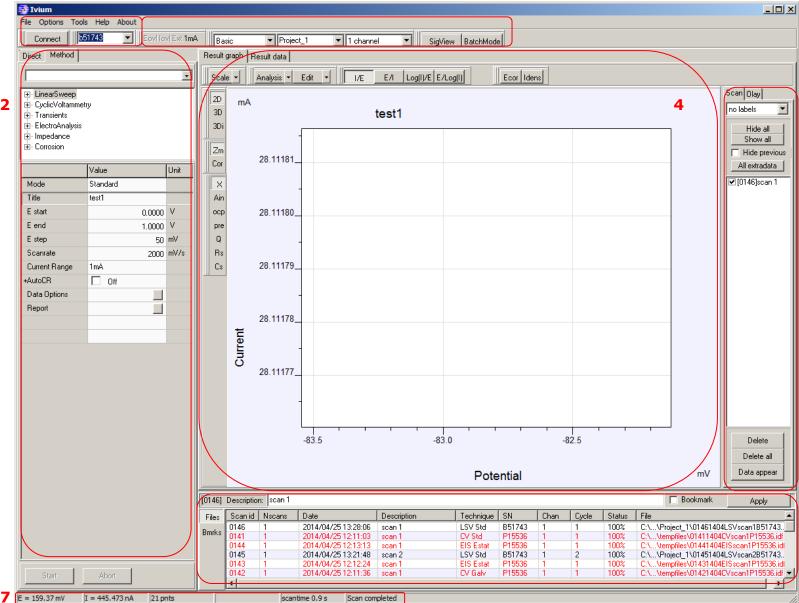

3. Advanced parameters

• In "Method" mode, by default the basic parameters are • The "Project window" shows the active Project directory shown. Selecting "Advanced" from the drop down menu in which data files are automatically stored. Select your will add a number of advanced method parameters. Project directory from the drop down menu. When typing a new Project name in the window, it will automatically be created when the next data file is stored.

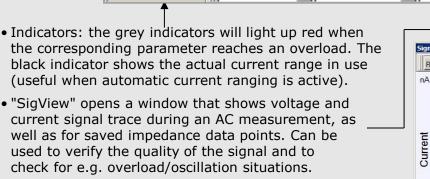

This can be used for automating and sequencing measurements, scheduling events, etc. For instructions, refer to the Technote "T1 Batch mode..." in your "IviumStat\Manuals and Release Notes" directory.

▼ 1 channel SigView BatchMode 0.2 -0.0 -0.2 -04 0.0 0.2 0.4 time Freq. = 1414.18 Hz Zabs = 3.389E+07 Ohm Phase = -86.32 deg status :0K(4:2:1)

Result graph | Result data | E scan |

0.05_

0.00

-0.05

Scale - Analysis - Edit - ZJPhi Rs, Cs Z', Z'' Y', Y'' Z,f

Result

5. Legend panel

Three tabs:

1) "Scan" shows the list of scans that have just been measured or loaded from memory. "Hide all"/"Show all" will respectively hide and show all scans in the list. In the list of scans, checking/unchecking the box next to the scan will show/hide individual scans. Note that the scan that is highlighted blue is the scan in active memory. This is the one that is evaluated in "Analysis", or has the numerical data shown in "Result data", etc. Clicking on a different scan will select that one (highlight it), making it available for analysis, saving, etc.

Checking the box before "Hide previous" will hide all previous scans, except the latest (useful when a large number of scans is listed). "All extra data" will show the extra data, for example OCP measurement, of all

scans in the list. "Delete" will remove the selected scan from the view screen only*.

"Delete all" will remove all scans from the view screen only*. *Data is always automatically stored in a datafile and cannot

be permanently deleted from the Legend panel "Data appear" will allow user to change the appearance of the data: lines,


colours, symbols. 2) "Olay" will allow the user to load data from file and overlay it so that it can be

compared to the scan data that was just recorded. Several scans can be overlayed

at the same time. All the same options apply as in the scan-tab. Overlayed data is not available for analysis. 3) The "Chan" tab shows when analog inputs are sampled (advanced method

parameter), it allows the user to show/hide analog channels in the second graph.

- A scan in the list can be selected (highlighted in blue), but not be visible in the graph, and vice versa.
- When different electrochemical methods are used, the data cannot always be represented in the same graph (i.e. CV and EIS data). For this reason, the format of the most recent scan in the Legend panel is automatically displayed and only compliant scans are shown. Non-compliant scans are hidden.
- Running a CV-scan will automatically clear the graph of all present scans when the next experiment is started; other methods will add the new scan to the Legend

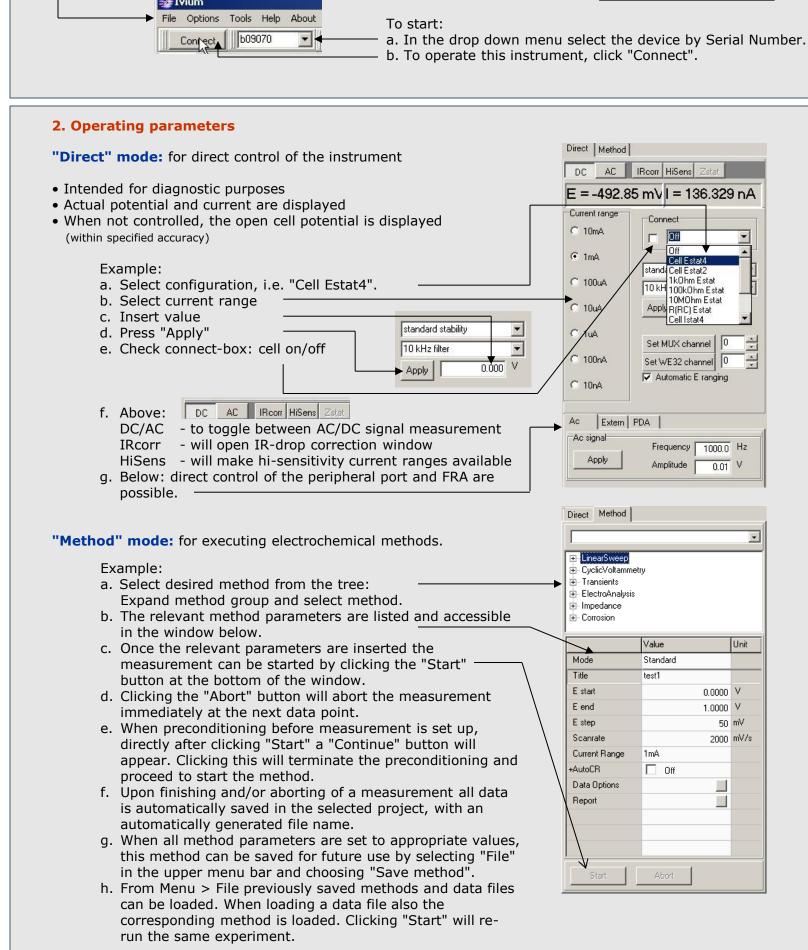

1) Install IviumSoft from CD:



IviumSoft will be placed on your desktop.




T2 Ivium Driver Installation.pdf

1. Device & software control

The menu bar:

C/ Software – user interface 1. Device & software control **Operating parameters**

4. Measurement results 5. Legend panel 6. Data file history list **Status bar**

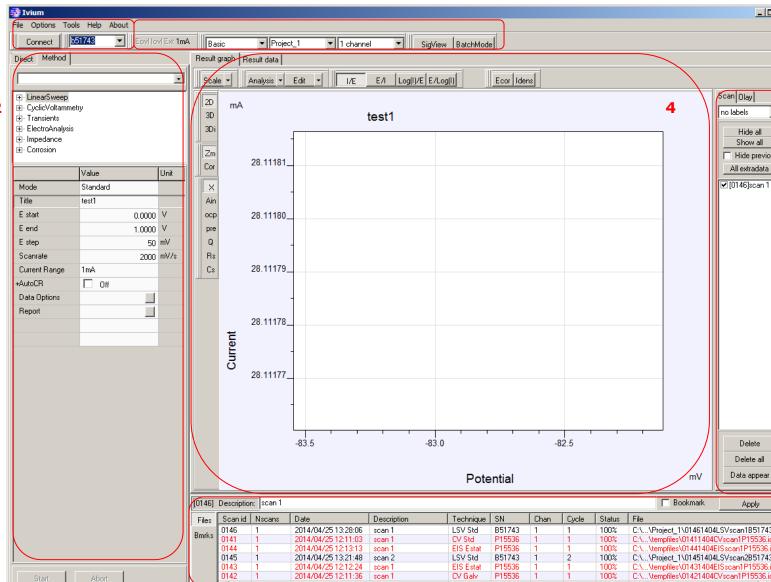

Eovillovi Ext 1mA ▼ Project_1 Basic • Indicators: the grey indicators will light up red when

• "BatchMode" opens the batch programming window.

4. Measurement results • When a measurement is started the data is shown real time in the result graph. A right mouse click on the graph will give the user a number of options, including changing the graph colour and background, as well as copying the graph to clipboard. The buttons to the left allow additional data to be displayed. Data is by default displayed "2D"; "3D" and "3Di" display may also be chosen. Clicking the "Cor" button after the measurement has finished, will start the correction mode, allowing the user to change individual data-points. "X" shows the primary data, "Ain, ocp, pre, Q, Rs, Cs" will all open a second graph that shows the corresponding data when available: optional analog inputs, ocpmeasurement before scan, pre-treatment data, etc.

- The "Scale" button allows the scale of the graph to be adjusted via the drop down menu; clicking on the button itself will auto-scale the graph.
- "Analysis" will make a number of analysis methods available to the data in active memory.
- "Edit" will allow smoothing of the scan data.
- The buttons at the top of the graph activate different representations of (impedance) measurement data.
- The tabs "Result graph" and "Result data" above the graph enable switching between graphical and numerical representation of the scan data. "E scan" will show the data for analysis when an impedance-scan method has been carried out.

6. Data file history list

All Ivium experimental data is automatically saved upon completion or user-abort of a measurement. A data file is created with an automatically generated file name, including a unique scan id and instrument S/N. After the data is saved to file in the active project, it will be added to the history list of most recent data files at the bottom of

> After an experiment is completed and the datafile is stored, the user can enter a description and click "Apply" to store the description in the view of the history list.

innovative solutions for electrochemical research

Checking the bookmark box and pressing "Apply" will bookmark the selected datafile for future reference. Bookmarked files are underlined.

File Options Tools Help About

Device maintenance

Performance tes

Pulse Generator

CurrentInterrupt

Test performance: connect Testcell 1 to cellc...

Connect TestCell 1, and press <Run test> to start

0.0000 V

0.0000 V

1.00 mA

1.000 mA

100.0 µA

10.00 µA

1.000 µA

100.0 nA

10.00 nA

1.000 mA

0.0000 V

1000.0 Ohm

Abort

1.0000 µF

100.0 Ohm 90.0

E_offset

I_10mA

I_1mA

I_100uA

1_10uA

I_1uA

I_100nA

1_10nA

I_cable

E_cable

AC_test_R1

AC_test_R2

AC_test_C1

Run test

Close

E_applied

Value Unit min limit max limit status

0.90

0.997

98.0

9.97

0.980

99.0

9.50

0.990

980.0

0.9970

-0.0030 0.0030

0.9970 1.0030

1.10

1.003

102.0

10.03

1.020

101.0

10.50


1.010

1.0030

300.0

1020.0

0.8000 1.2000

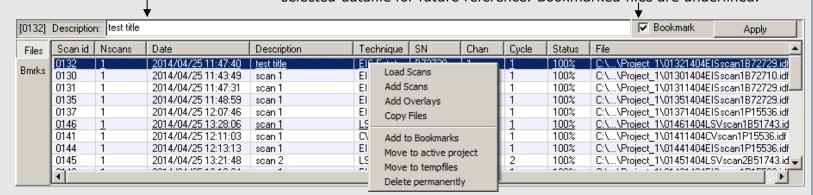

Double clicking on a file in the history list will load that file into the result window. Holding the [Shift] key will allow multiple data files to be selected simultaneously. A right mouse click will open a pop-up:

- Load Scans: will load selected scans into the result window.
- Add Scans: will add selected scans to the result window and the scans already there.
- Add Overlays: will load selected scans to the result window but into the Olay tab, in stead of the Scan tab.
- Copy files: will copy data file to clipboard, for example for easy pasting into an e-mail, or a file directory. • Add to Bookmarks: will add the selected scan to the bookmarks for future reference. Book marked scans can be identified by their underline. All bookmarked files in the history list can be shown by clicking [Bmrks] to the left
- Move to active project: will move the data file to the active project (see 3. Advanced parameters).
- Move to tempfiles: will move the data file to the tempfile directory. Data files in the "tempfile" project are displayed in red.
- Delete permanently: will delete the data file permanently.

Data Storage:

Data files are stored in a Library\project structure. In the menu "File>Data Explorer" data files in all libraries and projects can be explored, with a list of method parameters and result graph preview. In the menu "Options>Datahandling Options" the data storing options can be specified, such as: timed intermediate saving, automatic storage of a file copy to another designation (i.e. network drive), temp file properties, etc.

7. Status bar

The Status bar is located at the bottom of the user interface. Depending on the status of the instrument (idle, running a method, etc.) and/or the method that is running, a series of parameters is shown.

E = -0.38 mV I = -498.070 nA 11 pnts cycle 1 scantime 2.5 s Scan completed

In the case of a CV scan (see figure above) it shows the real time E, I, number of acquired data points so far, cycle number, elapsed runtime of the scan and the notification that a measurement is in progress (or completed). When, for example, an OCP measurement is included in the method, the measured final reference OCP value is also shown in the status bar. During an impedance measurement the real time E and I, the current frequency of the datapoint measured and the percentage of progress are displayed.

D/Special operations

From the top menu "Tools" will open a list of special operations: -

"Performance test" (only available when an instrument is connected)

"Performance test" will open a pop-up window that allows a self-diagnostic test of the instrument. This test is intended for occasional use, for example to determine if the calibration is still correct or whether the instrument is in proper working order. At a service request you can be asked to run this test by an Ivium service engineer to determine if the problem is hardware related. The performance test can only be executed when the instrument is in the "standard" configuration (= no extended range or other options activated).


To run the test, connect the cell cable to Testcell1, uncheck all options in the menu "Options>Options", and run the test. All status reports should read "Pass". If one, or more, reads "fail" after repetitive runs, the instrument may need calibration. This cannot be done by the user, you should contact your supplier.

The test report is stored in your\Iviumstat directory as a text

file with the file name: [Serial number].ipt

"Pulse generator"


"Pulse generator" is a direct function that will apply a pulse according to the selected parameters:

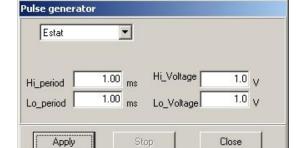

By selection, a potentiostatic or galvanostatic pulse can be generated with user selectable height and duration. The pulse generator is intended for preparation/conditioning of a sample. Data is not measured or stored.

"Current Interrupt"

"Current Interrupt" will open a window that allows use of the Ivium Current Interrupt module (CIM). The CIM can be purchased separately. The current interrupt technique can be used to determine the ohmic drop of an electrochemical system. This value can then be used for ohmic drop compensation during various electrochemical methods.

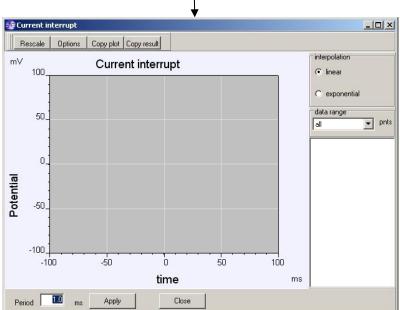

"Define channels"

"Define channels" is only available for the Ivium-n-Stat: it will open a pop-up window that allows the user to assign the channel serial number to a channel number for easier designation and reference. Details on how to do this can be found in the IviumSoft Help 'Getting started>Ivium-n-Stat'.

E/ Software & Firmware upgrade

IviumSoft upgrade

IviumSoft is frequently updated with new features, often at a customer's request. When a new version of IviumSoft is available, it is published on our website at: www.ivium.nl/Support. Here the new version can be downloaded as .zip file. The .zip file contains the IviumSoft.exe and whatever auxiliary files necessary. Note that this does NOT contain a full installation package, so using this file for a new installation is not possible.

To upgrade the IviumSoft, download the latest version of the software from www.ivium.nl/Support. Open the .zip file and copy all files to clipboard. Browse to your IviumStat file folder: ..\IviumStat, and paste (replace the files already there with the new ones).


Now your IviumSoft has been upgraded. After this proceed to upgrade the firmware of your instrument.

Firmware upgrade

Ivium instruments are equipped with an internal micro PC that operates the electronics, and is used to store data during a HiSpeed experiment. The firmware that runs on the micro PC is specific to the version of IviumSoft. For correct communication (and operation) these versions need to match, that is why IviumSoft has the correct version of firmware embedded. To upgrade the firmware of your instrument:

1. Connect to your instrument in the IviumSoft and go to "Tools>Device maintenance"

2. In the pop-up window click on "Firmware upgrade"

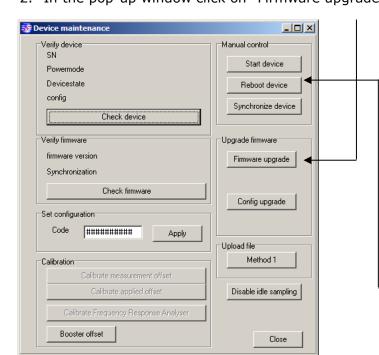

3. In the second pop-up click "Start", let the file upload, and when finished (successfully uploaded)

Connected

Direct Metho

⊕ LinearSwe

File Options Tools Help About

Device maintenance

Pulse Generator

CurrentInterrupt



- 4. After this, click "Reboot device".
- 5. Then close IviumSoft and restart your instrument.

For more instructions refer to Technote T4: "Software and Firmware Upgrade".

F/ Manual in Help-file

In addition to the brief overview of the IviumSoft user interface that is given in this Quick guide, in-depth information is given in the digital help file/manual that is included in your IviumSoft. There you will find instructions on the installation of potentiostats and modules, their instrument specifications, and how to operate them. In addition detailed information on how to use IviumSoft and all its functions regarding operating all electrochemical techniques and measurement methods, data representation and evaluation, etc., is described.

To access the digital help file/manual open your IviumSoft and press "F1". The help file is partly content sensitive so indicating with your cursor an area or subject in the IviumSoft user interface and subsequently pressing "F1" will open the helpfile on the relevant subject.

In the help file/manual the subjects are structured in main chapters according to:

Electrical Compliance Declaration of CE conformity of Ivium instruments.

Introduction Brief description of Ivium, our instruments and the working principle of IviumSoft.

🦫 IviumSoft Detailed description of IviumSoft: how to operate the software, instructions on the user interface, data representation and options, how to select and set instrument options and operations.

Getting started Instructions for first time installation: IviumSoft and Ivium equipment, instrument drivers and configuring your Windows system, product-specific information for all Ivium potentiostats and modules, how to connect the electrodes and internal

Description of how the user can control Ivium potentiostats through IviumSoft: Instrument control directly for diagnostic purposes or via user selectable electrochemical methods; how to program a sequence of events in the Batch programming mode; special

instrument functions; sequence of execution of the various commands by the potentiostat.

Electrochemical

integrated dummy cells.

Brief description of all available electrochemical techniques (for detailed descriptions techniques the electrochemical textbooks should be consulted), grouped by: Linear sweep, Cyclic voltammetry, Transients (CA, CP, ECN, MixedMode: versatile combination of transients and sweeps), Electroanalysis, Impedance, Corrosion.

Measurement Results How to access and represent the measured data (graphic and numeric), as well as the various available data evaluation techniques and possibilities.

Data files How and where data files are stored and accessed; the structure of projects and libraries; the handling of very large data files and automatic saving of data.

Special functions Special operations for (instrument) diagnostics and maintenance, preparatory

techniques and use of the peripheral port signals and functions.

Software upgrade Instructions for IviumSoft upgrade.

Control via Description on the software development driver and dll for control via other software other software such as Labview, C++, VB, etc.

Trouble shooting How to solve some of the most frequently encountered issues and how to restore your potentiostat.

Instrument Specifications of all Ivium instruments, modules, accessories and their compatibility.

Method parameters Detailed description of all method parameters.

specifications

It may be that when you open the helpfile the content is not visible, or you see a cryptic browser message like "Navigation to the webpage was cancelled" or "Action cancelled". This is caused by Windows and can be fixed simply by changing a setting: In Windows explorer navigate to your IviumStat directory where the help file is located. Right-mouse-click on the 'IviumSoft.chm' file and choose 'Properties' at the bottom of the list. This opens a pop-up. In the bottom panel of the pop-up click "Unblock" and OK. Close the pop-up window. Now the content in the IviumSoft help will be available.

The help file/manual covers current IviumSoft and instruments and operations. If you have specific operating or instrument specification questions regarding older versions of Ivium instruments, please contact Ivium.

A (printable) PDF-version of the digital IviumSoft Help is available in the 'Manuals and Release Notes' directory in the 'IviumStat' installation directory on your PC.

